

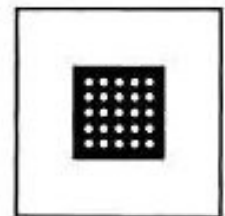
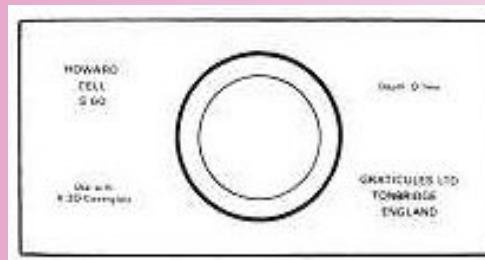
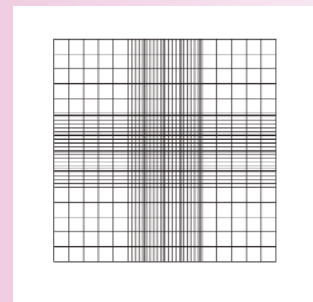
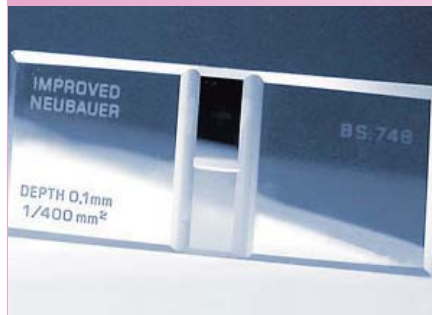


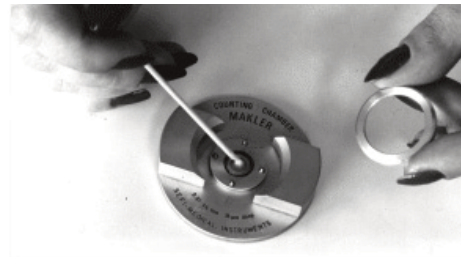

COUNTING CHAMBERS COUNTING CHAMBERS COUNTING CHAMBERS COUNTING CHAMBERS COUNTING CHAMBERS COUNTING CHAMBERS

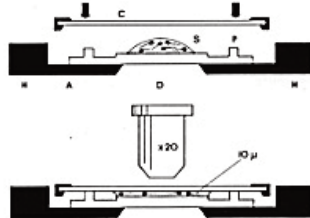

COUNTING CHAMBERS

As well as an extensive selection of specialist chambers for Haemocytometry, Parasitology and Cytology. The range includes:-

Pyser-SGI Counting Chambers.

The Pyser-SGI grid-marked Sedgewick Rafter counting chamber. Available in both plastic and durable glass versions, excellent for determining counts in known volumes of sample even using zoom stereo microscopes.

The Pyser-SGI Howard Cell for spore counting in tomato and fruit juice.

The MAKLER Semen Counting Chamber.



- Sedgewick Rafter Cell
- Howard Cell
- Makler Chamber
- Specialist Chambers

Pages 3-5

Pages 6-7

Pages 8

Pages 9-17

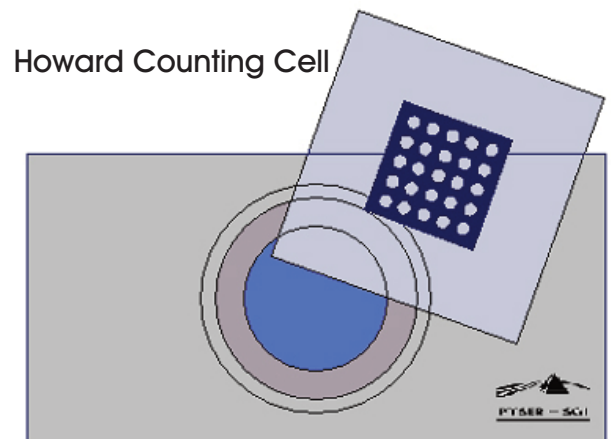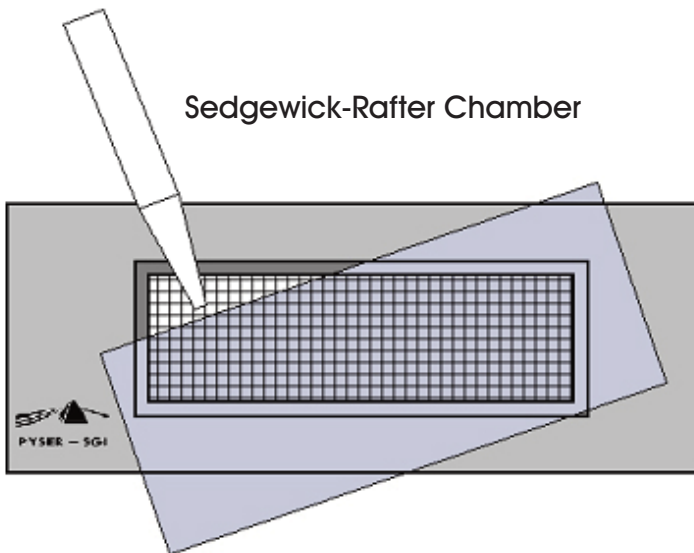

SEDGEWICK RAFTER CELL

The Pyser-SGI S50 & S52 Sedgewick Rafter counting cells are used widely in water analysis, culture inspection and for other applications where particles per unit volume in fluid must be determined.

Pyser-SGI's Sedgewick-Rafters are available in glass and plastic versions. The glass cell is designed for repeated cleaning and re-use, whereas the plastic version is designed for either a single or finite use only. Each chamber is supplied with one type S51 cover glass.

The Sedgewick-Rafter Chambers consists of a 2mm thick clear base slide (76 x 40 mm), onto which has been built a rectangular chamber. On to the top of this chamber is placed a cover slip.

This chamber is 50mm long x 20mm wide and 1mm deep and its base is marked with a grid of one thousand 1mm squares. When filled with liquid and with a cover glass placed over the chamber, a 1ml volume of liquid is trapped.

When used under a low magnification light microscope, each of the grid squares equates to 1microlitre (μ l) of liquid.

S50 - Plastic Sedgewick Rafter Counting Chamber

Pattern	Description	Order code
S50	Sedgewick Rafter Counting Chamber [plastic] Includes 1 cover glass	02C00415



Key Features

Plastic - economically priced cell


Description

An economically priced cell, for one time (or possibly several time) use and for educational/training exercises.

Generally we do not recommended the S50 plastic version for the professional user. The softness of the plastic means it is easily scratched and otherwise damaged. However most laboratories who do use the plastic version see it more as a "throw away" or consumable, which may best suit certain applications.

S52 - Glass Sedgewick Rafter Counting Chamber

Pattern	Description	Order code
S52	Sedgewick Rafter Counting Chamber [Glass] Includes 1 cover glass	02B00417

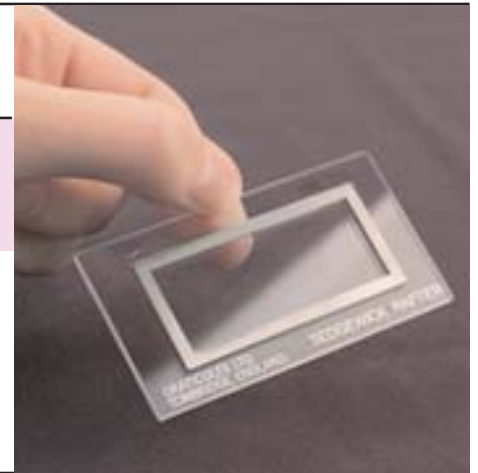

Key Features


Glass - Professional Cell

Description

A serious reusable tool for the professional user. The cell is made of high quality optical glass with a chromium surface image. It is intended for continued professional use and whenever using phase contrast.

S51 - Spare Coverglass for Pyser-SGI Sedgewick Rafter Counting Chamber

Pattern	Description	Order code
S51	Cover glass for use with both Glass & Plastic Cells	02C00416

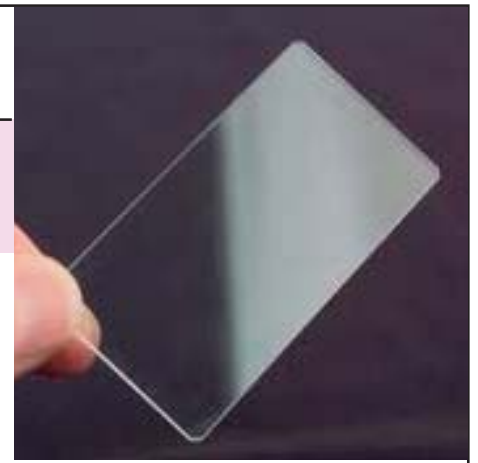

Key Features

For use with both Glass & Plastic Cells

Description

Both Pyser-SGI Plastic and Glass Sedgewick Rafter Cells are supplied with a cover glass. Additional spare cover glasses are available to order from stock, as this S51

Sedgewick-Rafter Filling Method & Usage

The cell can be used with either living or preserved material. To fill the cell, place the cover glass across the chamber top (Fig A). This allows the air bubbles to escape during the filling procedure. The sample is then taken into a 1ml wide mouthed pipette and then carefully transferred to the chamber. Do not overfill the chamber, because the volume of the sample in the chamber must be known exactly and the cover glass must not float free, but held onto the cell walls by surface tension. During counting, water may evaporate from the chamber. To prevent gas bubble formation, a small drop of distilled water may be placed on the slide outside the cell, just touching the cell wall and cover glass. Before the cell count is made the Sedgewick-Rafter chamber should be allowed to stand for at least 15 minutes to allow algae, or other particles, to settle to the bottom.

The Grid Pattern in the base of the chamber assists the counting and calculation process, by clearly defining a known sample volume in $1\mu\text{l}$ blocks

Counting in strips is easier, no need to use reticle grids or know the precise area of your field of view.

This makes Pyser SGI's S50 & S52 Gridded Sedgewick-Rafter Cell perfect for use with a Stereo Zoom Microscope without the need to carefully set the precise zoom/magnification before counting the sample.

For precise sample preparation and calculation methods, you should follow your own internal or published procedures.

Filling a Sedgewick-Rafter Chamber

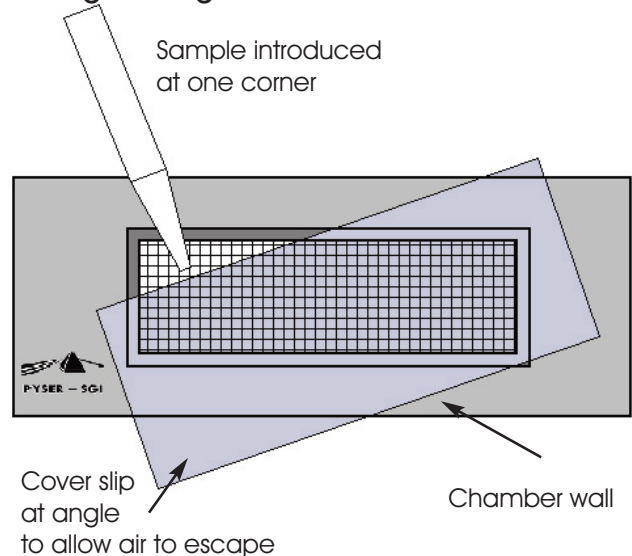


Fig A

Cleaning Method

To clean the counting chamber: After completing the count, remove the cover glass and clean the counting chamber with water or a mild cleaning solution (10% solution of bleach). Dry the counting chamber with a soft cloth or wipe, or rinse with acetone.

HOWARD MOULD COUNTING CELL

The Pyser-SGI S60 Howard Cell and associated K20 Cover Glass, is used world wide for mould counting in food quality control applications, such as tomato products and other fruit based preparations.

Pyser-SGI's improved Howard Cell is based on the method originally developed by B.J. Howard in 1911, primarily for the purpose of monitoring tomato products, using a microscope.

The special S60 Howard Cell chamber slide in conjunction with the K20 Cover Glass is designed to count mould mycelia.

The S60 Howard Cell is a glass slide 76mm x 35mm with a central circular island and is used for counting mould fibres and spores in fruit juices especially from tomatoes. With the K20 cover glass in place a 0.1mm thickness of liquid is contained over the central island.



The cover glass has 25 calibrated fields of 1.382mm diameter through which to view the particles.

This cover glass removes the necessity of precise adjustment of the microscope magnification and calibration of a special eyepiece reticle in the original Howard Method, making it suitable for use with a modern zoom stereo microscope as well as a conventional compound microscope.

Note: The complete system requires both the cell itself and cover glass. The cell does not come with a cover glass and one must be ordered separately. We recommend ordering one or two extra cover glasses, this way there is always going to be one in reserve.

S60 - Howard Cell for Fruit Juices

Pattern	Description	Order code
S60	Howard Cell for Fruit Juices	02C00419
K20 Cover Slip Required for use		

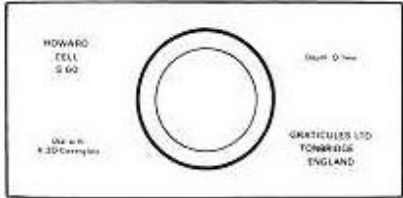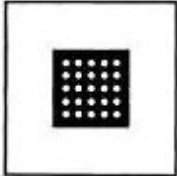Key Features


Special Chamber for use with K20 windowed cover slip to count mould fibres and spores

Description

The S60 Howard Cell is a glass slide 76mm x 35mm with a central circular island and is used for counting mould fibres and spores in fruit juices especially from tomatoes. With the K20 cover glass in place a 0.1mm thickness of liquid is contained over the central island.

K20 - Cover Glass for Pyser-SGI Howard Cell

Pattern	Description	Order code
K20	Coverglass for Howard Cell	02C00420

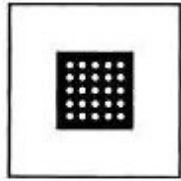

Key Features

For use with S60 Cell to hold a 0.1mm thickness of liquid for analysis and present known sample volumes for counting

Description

When used with the S60 Howard Cell, the K20 cover glass holds in place a 0.1mm thickness of liquid is over the central island for analysis.

The cover glass has 25 calibrated fields of 1.382mm diameter through which to view the particles.

This cover glass removes the necessity of precise adjustment of the microscope magnification and calibration of a special eyepiece reticle in the original Howard Method, making it suitable for use with a modern zoom stereo microscope as well as a conventional compound microscope.

How to Use a Howard Cell - General Procedure

The material to be examined should be a pulp. Mix a small quantity with water until the solids of the diluted pulp are between 8.37% and 9.37%. This corresponds to an Abbe refractometer reading at 20°C of 1.3460.

Spread a small drop of the well-mixed sample with the end of a glass rod over the counting chamber.

Place the cover glass on to the counting chamber and carefully press down the shoulder of the chamber until Newton's rings are visible.

Prepared samples containing air bubbles beneath the cover glass or an over-full moat should be discarded.

If using a compound microscope, examine using the X10 eyepiece and the X10 objective.


Systematically examine all 25 fields and note those with a presence or absence of mould filaments (hyphae).

A field is regarded as positive if the aggregate length of not more than three filaments present exceed one sixth of the diameter of the field.

This is a general description of how a Howard Cell is used. The results are interpreted as a percentage of positive fields observed in all the fields examined.

Precise interpretation of the results is made by a statistical analysis of the sample and should be carried out in accordance with your own internal or published procedures.

Howard Mould Counting Cell & Cover Glass

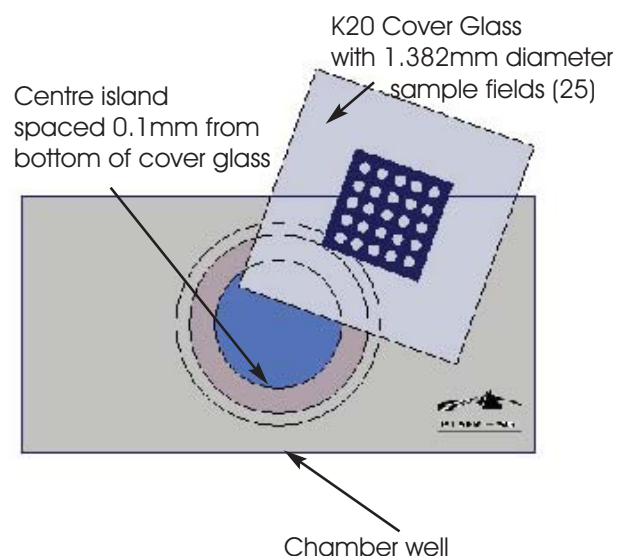

MAKLER CHAMBER



The Makler counting chamber is only 10µm deep, which is one tenth the depth of other Haemocytometers, making it the shallowest of known chambers.

This quality chamber is constructed from two pieces of optically flat glass: the first is the chamber, the second the cover glass which has a fine grid pattern of 1mm squares and a centre area further subdivided into 0.1mm squares.

Four quartz pins of precise height hold the cover glass to give an exact trapped specimen depth of 10 µm

Makler Chamber

Pattern	Description	Order code
Makler	0.01mm deep Semen Counting Chamber	02C00645

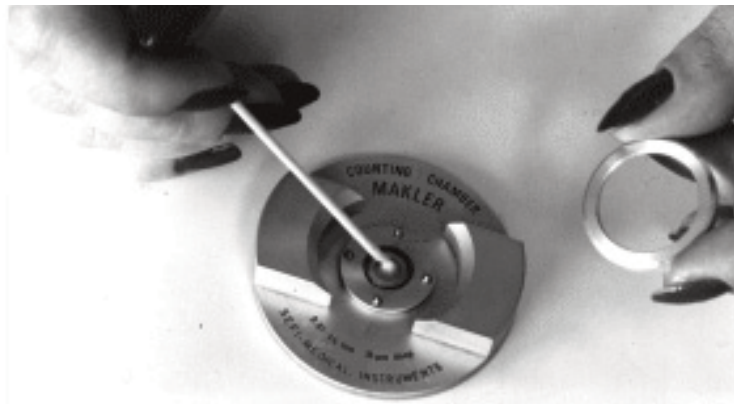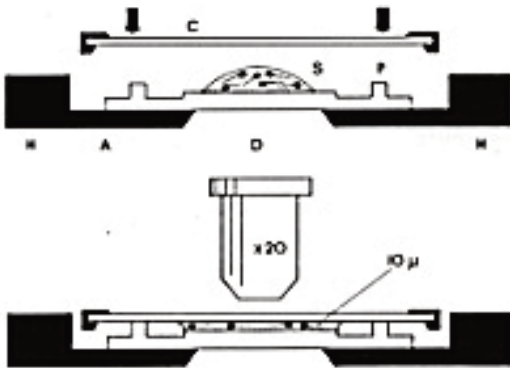

SPECIALIST CHAMBERS

This range of quality counting chambers are available in a wide range of types for Haemocytometry, Parasitology, Cytology and other applications

Chambers with inside markings have rulings applied using engraving techniques to give a permanent image, clearly visible through the glass without sacrificing light transmission.


Chambers with ruled cover glass have robust image applied using chrome deposition.

Haemocytometry		page
Agasse-Lefont	Single Cell	12
Burker	Double Cell	12
Malassez	Double Cell	10
Malassez	Single Cell	10
Neubauer	Double Cell	11
Neubauer	Single Cell	11
Neubauer, Improved	Double Cell	12
Thoma	Double Cell	10
Thoma	Single Cell	11
Turk	Double Cell	13
Parasitology, Cytology & Others		page
Agasse-Lefont B	Single Cell	16
Evaluation	Single Cell	15
Fertility	Single Cell	16
Helber	Single Cell	15
Lemaur	Single Cell	14
McMaster, Improved	Double Cell	15
Nageotte	Double Cell	14
Nageotte	Single Cell	14
Chamber Cover Glasses		page
McMaster, Old	Cover Glass	17
22x22mm x0.45mm	Cover Glass	17
30x28mm x 0.45mm	Cover Glass	17

Haemocytometry Counting Chambers

Malassez Double Cell

Pattern	Description	Order code
Malassez	Double Cell 0.2mm	02C00600

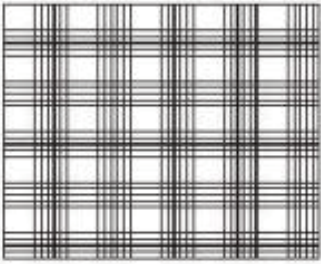


Key Features
For Haemoglobin and Lucocyte counts

Description
Standard French Ruling

Malassez Single Cell

Pattern	Description	Order code
Malassez	Single Cell 0.2mm	02C00601

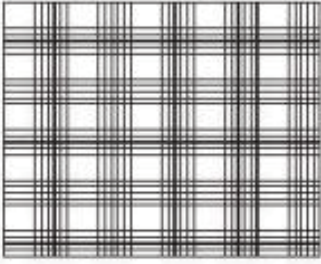
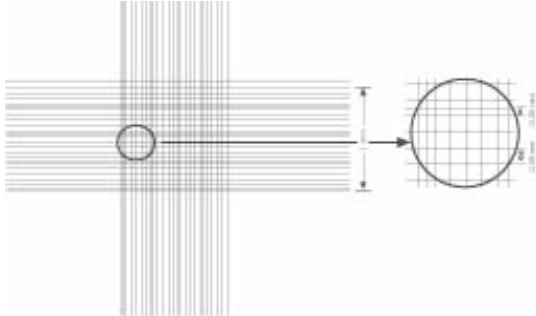


Key Features
For Haemoglobin and Lucocyte counts

Description
Standard French Ruling

Thoma Double Cell

Pattern	Description	Order code
Thoma	Double Cell 0.1mm	02C00605

Key Features
For Haemoglobin and Lucocyte counts

Description
Standard German Ruling


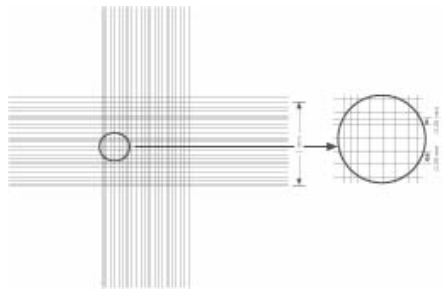
Haemocytometry Counting Chambers

Thoma Single Cell

Pattern	Description	Order code
Thoma	Single Cell 0.1mm	02C00606

Key Features
For Haemoglobin and Lucocyte counts

Description
Standard German Ruling


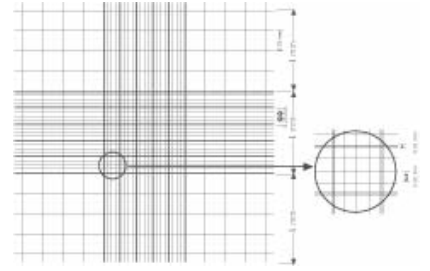



Neubauer Double Cell

Pattern	Description	Order code
Neubauer	Double Cell 0.1mm	02C00610

Key Features
For Haemoglobin and Lucocyte counts

Description
Modified Thoma Ruling


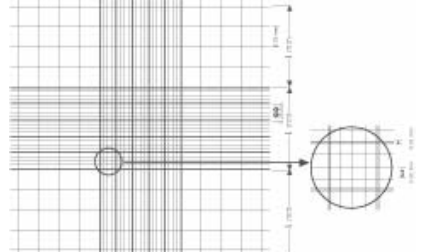



Neubauer Single Cell

Pattern	Description	Order code
Neubauer	Single Cell 0.1mm	02C00611

Key Features
For Haemoglobin and Lucocyte counts

Description
Modified Thoma Ruling

Haemocytometry Counting Chambers

Improved Neubauer Double Cell

Pattern	Description	Order code
Improved Neubauer	Double Cell 0.1mm	02C00616




Key Features

For Haemoglobin and Lucocyte counts


Description

Ruling allows full use of central squares

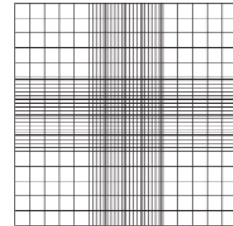

Burker

Pattern	Description	Order code
Burker	Double Cell 0.1mm	02C00626

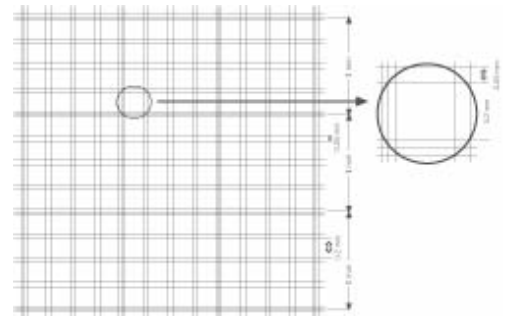

Key Features


For Haemoglobin and Lucocyte counts

Description

For counts in bands

Agasse-Lefont R

Pattern	Description	Order code
Agasse-Lefont R	Single Cell 0.001mm	02C00670

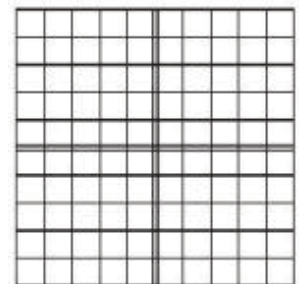

Key Features

For Haemoglobin and Lucocyte counts

Description

Completely Decimal



Haemocytometry Counting Chambers

Turk Double Cell

Pattern	Description	Order code
Turk	Double Cell 0.1mm	02C00621

Key Features
For Haemoglobin and Lucocyte counts

Description
Combination of Burker and Thoma Rulings

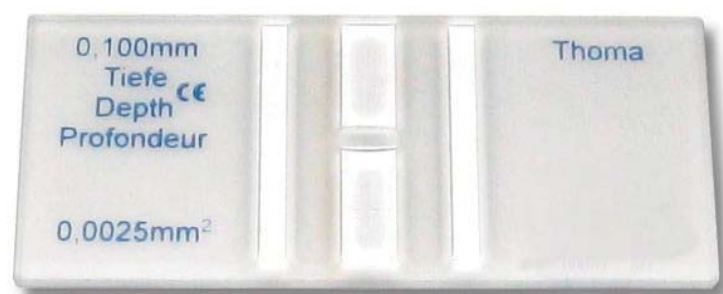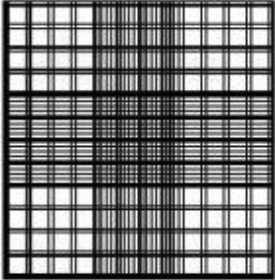

Typical Counting Chamber


Parasitology, Cytology & Other Chambers

Nageotte Double Cell

Pattern	Description	Order code
Nageotte	Double Cell 0.5mm	02C00630

Key Features
For Urinary cytology, cephalo-rachidien fluid analysis

Description
For liquids in poor content (French)

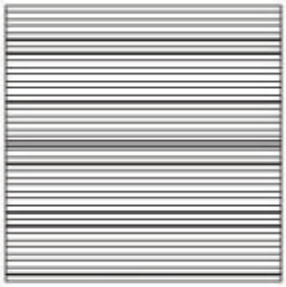


Nageotte Single Cell

Pattern	Description	Order code
Nageotte	Single Cell 0.5mm	02C00631

Key Features
For Urinary cytology, cephalo-rachidien fluid analysis

Description
For liquids in poor content (French)

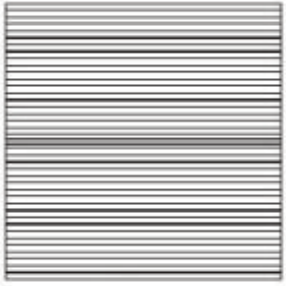


Lemaur Single Cell

Pattern	Description	Order code
Lemaur	Single Cell 0.4mm	02C00635

Key Features
For Urinary cytology, cephalo-rachidien fluid analysis

Description
Improved Nageotte with smaller volume



Parasitology, Cytology & Other Chambers

Improved McMaster Double Cell

Pattern	Description	Order code
McMaster	Double Cell 1.5mm	02C00650




Key Features
Parasitology

Description
Worm egg counting

Chamber is ruled with 2 x 1cm squares, each with 10 equal distance vertical lines forming 10 mm squares with 10 equal vertical columns.

Helber Single Round Cell

Pattern	Description	Order code
Helber	Single Round Cell 0.02mm	02C00655

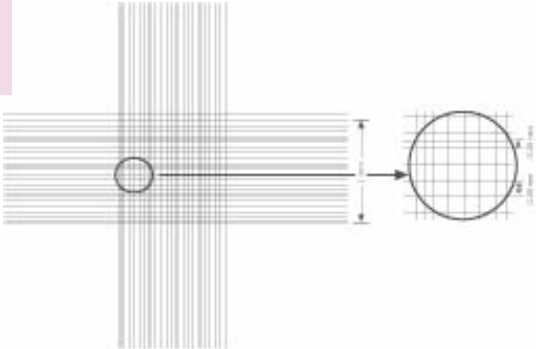

Key Features
Bacteria

Description
Slide 1mm thick with Thoma Ruling

Evaluation Single Round Cell

Pattern	Description	Order code
Evaluation	Single Round Cell 0.02mm	02C00660

Key Features
Bacteria

Description
Slide 1mm thick - no ruling


Parasitology, Cytology & Other Chambers

Semen Single Round Cell

Pattern	Description	Order code
Fertility	Semen Single Round Cell 0.02mm	02C00665

Key Features
Fertility

Description
For counting untreated semen. Ruling of 10 x 10 of 0.1mm squares

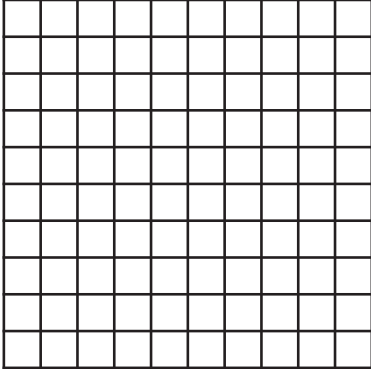


Agasse Lafont B Single Cell

Pattern	Description	Order code
Agasse-Lefont B	Single Cell 0.003mm	02C00675

Key Features
Other


Description
Completely decimal

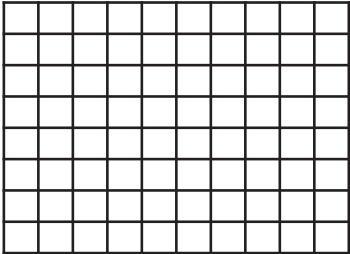

Chamber Cover Glasses

Cover Glass for old type McMaster

Pattern	Description	Order code
Old type McMaster	Cover Glass	02C00651




Key Features
Parasitology

Description
This cover slip is for the old McMaster chamber type see Improved McMaster for current product

22mm x 22mm Thickness 0.45mm

Pattern	Description	Order code
Pack of cover glasses	22mm x 22mm	02C00700

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


Key Features
Suitable for counting chambers size 22mm x 22mm or 22mm x 25mm

Description
Pack contains 5 pairs (10) cover glasses

30mm x 28mm Thickness 0.45mm

Pattern	Description	Order code
Pack of cover glasses	30mm x 28mm	02C00701

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

Key Features
Suitable for larger counting chambers

Description
Pack contains 5 pairs (10) cover glasses

Notes


Other products

Inspecta



Other Products in the Inspecta Range:

- Magnifiers with Reticles
- Thru Hole Scopes
- Microscope Components
- Calibration Products
- Portable Microscopes
- Microscope Systems
- Video Inspection
- Co-ordinate Measurement
- Depth Scopes
- Measuring Microscopes

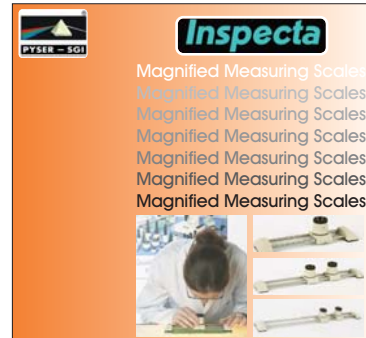

Other products



Please contact Pyser-SGI Limited



Pyser-SGI Limited reserves the right to amend these specifications.
 Visit our website on www.pyser-sgi.com for further information on the complete range of products available from Pyser-SGI Limited.



PYSER - SGI LIMITED
 Fircroft Way, Edenbridge, Kent,
 United Kingdom, TN8 6HA
 Telephone: +44 (0)1732 864111
 Facsimilie: +44 (0)1732 865544
 Email: sales@pyser-sgi.com
www.pyser-sgi.com
 July 2011 E & OE